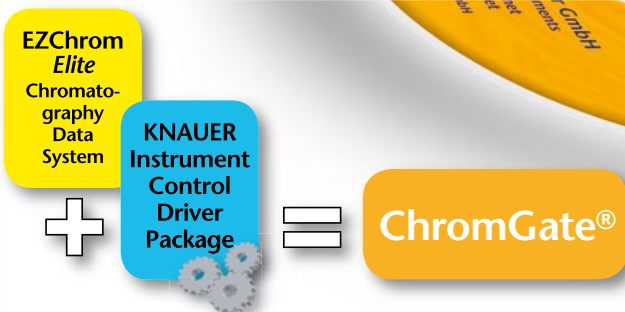
# Software ChromGate®

## HPLC software, not just for HPLC ...

ChromGate® is a full-featured scalable chromatography data system in compliance with FDA 21 CFR Part 11 for instrument control, data acquisition and report creation. It can be used to control a wide range of chromatography systems ranging from capillary electrophoresis to analytical and preparative HPLC, and even up to simulated moving bed systems. Chromatography tasks can be easily scaled-up while staying in **one single software environment**.

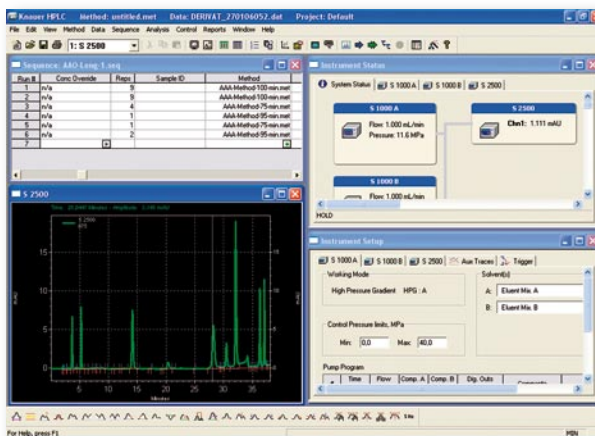
Benefits for the user are a unified user interface and simplified operation made possible by "wizard"-assisted programming. Lab managers benefit from reduced training expenses, a uniform system configuration and data reporting, as well as less maintenance.

ChromGate® software is based on the well-known EZChrom Elite kernel and the "Rapid-Control" open driver concept which makes it possible to incorporate and control non-KNAUER instruments into a ChromGate system. KNAUER instruments can also be integrated into other EZChrom Elite systems supporting Rapid-Control drivers. **More than 300** EZChrom Elite drivers for **instruments** from different manufacturers such as Agilent, Varian or Thermo Electron are currently available. With the ChromGate® Generic Driver Option, exclusively from KNAUER, virtually any instrument which communicates via RS-232 (ASCII) can also be integrated.

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
## The best solution for large and small laboratories

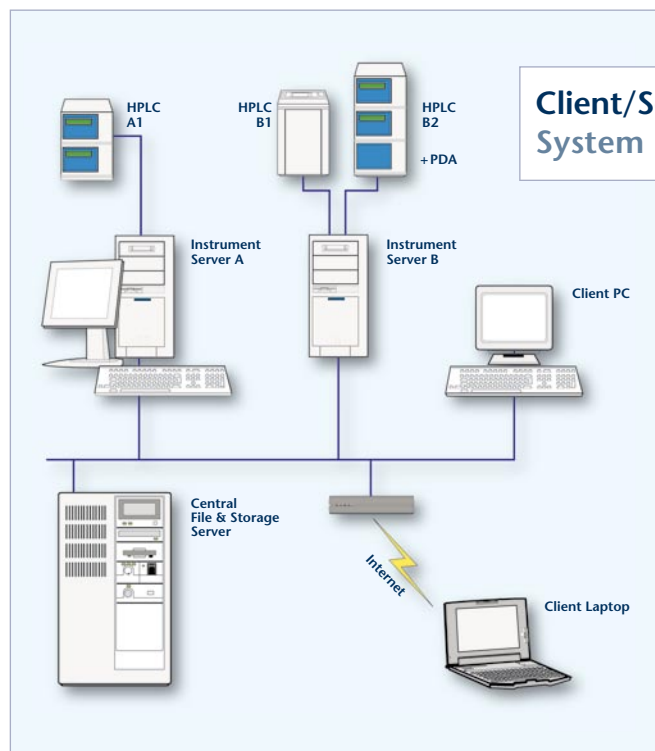Scalable from a single-user system to an enterprise wide Client/Server configuration, ChromGate® can deliver the performance required for large networked laboratories, yet is also compact enough to be used effectively in small laboratories.


The logical Client/Server architecture allows for ChromGate® to be accessed from every computer in the network. With ChromGate® it is possible to process data in the office as well as to control the HPLC system over the Instrument Server PC in the laboratory. Users and groups as well as access rights can be easily modified from the Windows™ domain controller.

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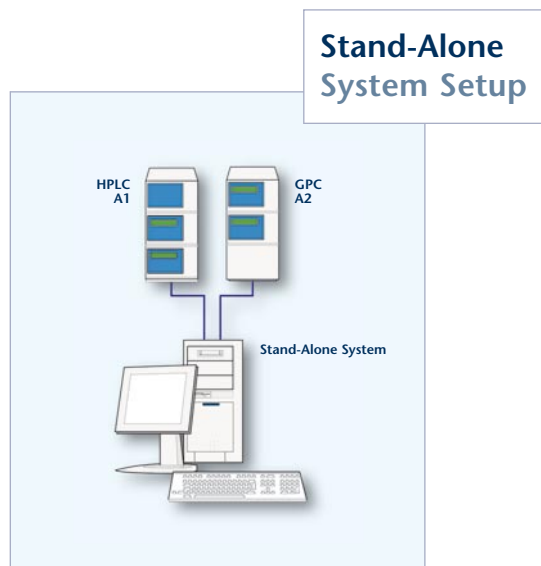### FDA 21 CFR Part 11

ChromGate® provides the highest level of data security available along with total GLP compliance. It fully implements the guidelines specified by FDA 21 CFR Part 11 for the storage and protection of electronic archived data and electronic signatures.

Other features include password-protected log in with definable access rights on the operating system level (Windows™ 2000/XP Pro), protection from inadvertent data deletion, as well as complete data traceability.

The method and the hardware configuration are saved with every data record so that all relevant information is accessible at any time. More features include an activity log and the Audit Trail Option for methods, sequences, and chromatograms.

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**Options**

Workstation | Enterprise | Email | General

General options

- Show all analysis results
- Automatically enable method audit trail for all projects
- Automatically enable password audit trail for all projects
- Activity log purge authorized only after archive
- Enable system activity log

Logging options

Security options

- SP: Expanded security
- Log out current user when idle for: [ ] min.
- Cancel an in-progress electronic signature when idle for: [2.0] min.

**Instrument Activity Log**

User	Logged	Source	User Name	App/ID
maine	18.12.2005 17:02	SERGEI_TEST	User Name logged in	
maine	18.12.2005 17:18:11	SERGEI_TEST	Run Queue - Complete Single Run - IL4	
maine	18.12.2005 17:13:43	SERGEI_TEST	Run Queue - Start Single Run - IL4	
maine	18.12.2005 17:13:42	SERGEI_TEST	Run Queue - Add Single Run - IL4	
maine	18.12.2005 17:13:06	SERGEI_TEST	Run Queue - Complete Single Run - IL4	
maine	18.12.2005 17:09:28	SERGEI_TEST	Run Queue - Start Single Run - IL4	
maine	18.12.2005 17:09:27	SERGEI_TEST	Run Queue - Add Single Run - IL4	
maine	18.12.2005 17:07:32	SERGEI_TEST	User Name logged in	
maine	18.12.2005 16:47:57	SERGEI_TEST	User Name logged out	
maine	18.12.2005 16:18:01	SERGEI_TEST	Direct Control: S 2000 Set HPLC Lang	
maine	18.12.2005 16:17:59	SERGEI_TEST	Direct Control: S 2000 Set D2 Lang	
maine	18.12.2005 16:22:49	SERGEI_TEST	Run Queue - Complete Single Run - IL4	
maine	18.12.2005 16:22:47	SERGEI_TEST	Run Queue - Add Single Run - IL4	
maine	18.12.2005 16:22:47	SERGEI_TEST	Run Queue - Start Single Run - IL4	
maine	18.12.2005 16:22:47	SERGEI_TEST	Run Queue - Add Single Run - IL4	

**Login**

User name: [Matthew]  
 Password: [ ]  
 Project: [Customize] [ ]

**Set User Privileges**

What privileges should this user have in each project?

User: [Testlaborant]

Selected projects: [Kundenkonto: JCP/MS]

Unassigned privileges:

- Method
- Save Method
- Advanced
- System Suitability
- Data
- Save Data
- Electronic Signature
- Sign Data Files

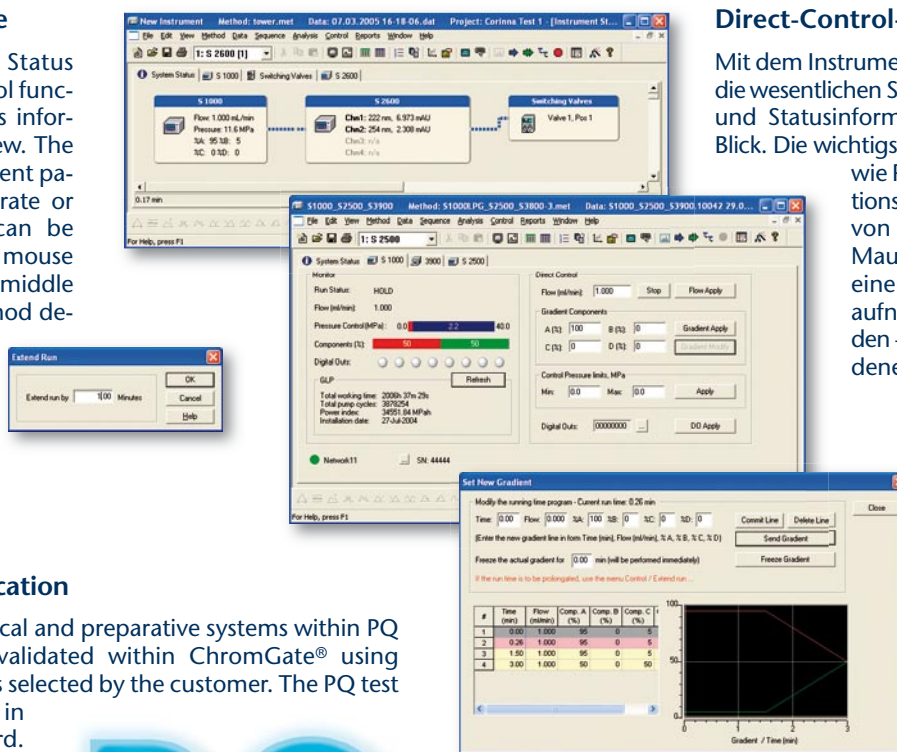
Assigned privileges:

- Open Method
- Instrument Setup
- Integration Events
- Peak Groups
- Custom Report
- Review Calibration
- Calibrate
- Data Files

## ChromGate® sets the standard for Chromatography Data Systems

### Direct-Control Mode

With the Instrument Status window, essential control functions and current status information are always in view. The most important instrument parameters such as flow rate or detector wavelength can be adjusted with a single mouse click, even while in the middle of a run – ideal for method development.

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### Performance Qualification

The operation of analytical and preparative systems within PQ specifications can be validated within ChromGate® using methods and conditions selected by the customer. The PQ test procedure is supported in the software as a wizard.

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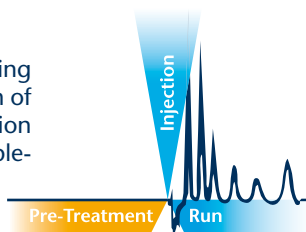
### Pre-Treatment Programming

The Pre-Treatment function allows for programming and control of pumps and valves prior to injection of the sample. With this function sample preparation or pre-run column conditioning steps can be implemented.

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
### Solvent Control

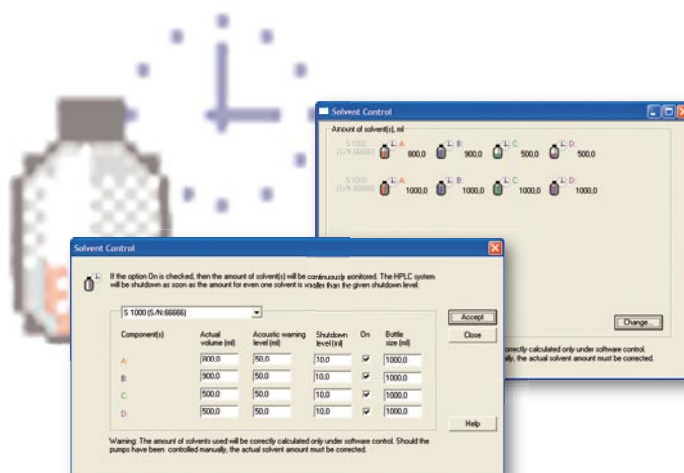Has your HPLC system ever run dry? ChromGate® does the thinking for you and calculates the expected eluent consumption with the Solvent Control feature for Smartline instruments. You can establish by what amount of solvent remainder you want to be warned and when the pumps should be shut off.

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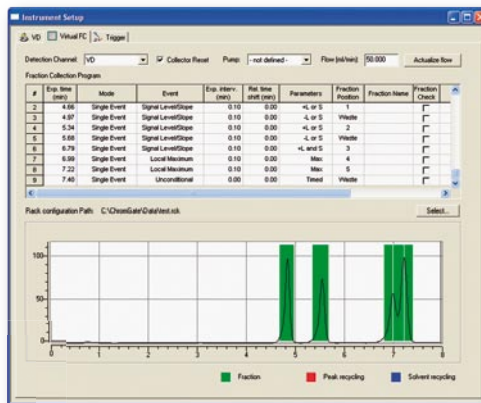
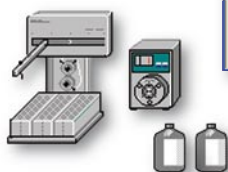
## ChromGate® Options for Support of Special Applications

### Preparative Chromatography

The Fraction Collector Control (FRC) option includes valuable tools for detector controlled fraction collection, solvent recycling and peak shaving.

Fractionation can be controlled by time (volume), level, slope, and spectra comparison with AND/OR linking of the switching criteria. Optimization of the fractionation is possible using only one master chromatogram, sparing additional need for sample and solvent. Some of the other remarkable features include an auto-recover option for interrupted runs and a direct control mode for interaction with running methods. The FRC Option transforms ChromGate® into a full-featured preparative chromatography software.

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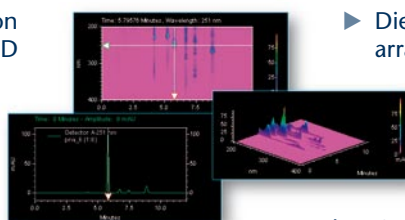
### Other ChromGate® Options

- ▶ With the **PDA** (photo diode array) option you can take full advantage of 2-D and 3-D chromatograms, peak purity analysis, or spectrum search in a spectra library.
- ▶ A **GPC/SEC Option** for molecular weight determination in size exclusion chromatography with various calibration methods.
- ▶ The **System Suitability Option** automates the calculation of system suitability parameters for system validation.

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Contact your local distributor ([www.knauer.net/worldwide](http://www.knauer.net/worldwide)) for more information on how ChromGate® can benefit your laboratory and request your free 7-day trial version.

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## Technical Data

### System

32-bit Chromatography Data Management System with Client/Server technology for instrument control, acquisition of data, data processing, and report creation under Microsoft Windows™ 2000/XP Pro.

### Expandability

**Stand-Alone:** One PC carries out all tasks from system control to data storage, to data analysis and reporting.

**Client/Server:** The HPLC system(s) is connected to an Instrument Server over which system control and data acquisition take place. Programming and data processing can also be carried out from any client PC in the network or even via the internet. Data can be stored on a central file server, making backups effortless. Automatic local buffering protects against data loss in case of IT network disruptions.

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### Supported instruments

More than 300 with available EZChrom compatible RapidControl drivers. With the flexible ChromGate® Generic Driver Option, the software can be adapted to even more instruments with RS-232/ASCII communication. (Please inquire.)

### Instrument Connection

One instrument (or up to 2 KnauerNet instruments) per RS-232 interface. PDA detectors either by Ethernet or PCI card. Non-KNAUER instruments directly or via KNAUER interface module.

### PC Hardware Recommendation

Min. Pentium IV 1500 MHz; 256 MB RAM (depending on operating system), additional 128 MB required per additional HPLC system or PDA detector; > 10 GB disk space.

### Data Import & Export

ASCII, ANDI standard, and OLE (object linking and embedding) for import into documents of other applications; ODBC for exchange with Windows™-based LIMS or databases.

### Graphics Capabilities

Multiple chromatogram view: zoom, print, cut, paste, copy. Mathematical operations such as align, stretch, move, or stacking of chromatograms. Modification of line color, type, and thickness. Full control of annotations for all peak parameters (e.g. angle, font style, size, and color).

### Integration

Real-time analysis of peaks, automatic or manual integration and baseline correction.

### Security and GLP

System integrity test of the software by validation tool. FDA 21 CFR Part 11 conformance, control of user privileges, various access levels, method, sequence, and data audit trails, electronic signature. Method parameters are saved in original data file and are protected against manipulation through checksums. Activity log files for instruments.

### Instrument Control

Instrument status display and Direct Control Mode are also possible while acquiring data, "wizards" simplify programming of methods and sequences. SmartSequence™ feature allows for automatic execution of pre-programmed functions (such as re-inject, stop, e-mail notification, etc).

### Calibration

Clearly-presented, editable calibration table. Interactive peak selection by mouse. Calibration curve fitting. Removal of outliers. Automatic peak naming and ordering.

### Report Function

Custom reports with an extensive set of options. Predefined templates.

### Other Features

Solvent level control for Smartline instruments. Performance Qualification wizard, pre-treatment.

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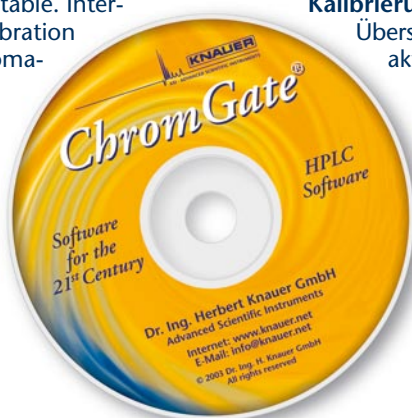
## Ordering Information

KNAUER offers brand-name PCs on which ChromGate® has been completely preinstalled and tested. Our PCs meet all ChromGate® requirements and are at the current state of the art when delivered.

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The CE (capillary electrophoresis) Option and the SMB (simulated moving bed) Control Option are available only when ordered with a KNAUER Capline System or KNAUER CSEP® SMB System, respectively.

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Please ask about our cross-upgrade discount!

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Dr. Ing. Herbert Knauer GmbH